




**FOR A SANITARY ENVIRONMENT  
IN ALL WORKPLACES**

# Lumitester™ Smart

Hygiene monitoring  
in three easy steps

Access from anywhere  
with cloud storage

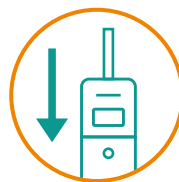

Visualize the data with App

With A3 Detection

**ATP+ADP+AMP**

**More Reliable and Sensitive | A3 Finds What Others Miss**

....

HyServe



# MAKE HYGIENE MONITORING A HABIT FOR EVERYONE

With the App, anyone can check cleanliness easily




## Measure | Easy-to-use, results within 10 seconds

STEP 1

### Swab


Swab the sample and insert the swab into main body

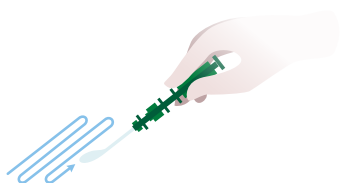

STEP 2

### Measure

Shake well and insert into the Lumitester

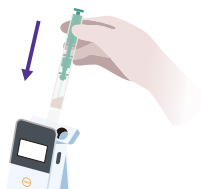

STEP 3

### Analyze

Measurements (RLU) displayed within 10 seconds



## Understand | App to continuously monitor multiple data

Recorded data can be displayed in time-series by registering with our specialized App. Inspection pass rates are automatically graphed and improvements can be visualized. Employees will be more conscious of hygiene, and will maintain a high standard of environmental sanitation, which will help build greater trust in your business



Display time-series data for each inspection



Visualize overall inspection scores through graphs



## Connect | Centralized multi-site data with cloud data management

Data will be stored in the cloud and can be accessed from anywhere in the world. This will allow fast detection and response to any problems that may arise.



### ■ Lumitester™ Smart Product code: 1902653

Measurement time	10 seconds
Data output	RLU (Relative Light Unit)
Power	2 alkaline or nickel hydride rechargeable batteries (AA)
Accessories	2 alkaline batteries (AA), Cleaning brush, USB cable, Strap, Quick Manual

※ Do not use this product for purposes other than hygiene monitoring.  
 ※ Lucipac A3 and Lumitester Smart should not be used for counting general living bacteria or detecting specific pathogens.

LuciPac™ A3 Surface	Product code: 1702671	100 swabs/kit
LuciPac™ A3 Water	Product code: 1702672	100 swabs/kit

### Storage Condition 2-8°C (Do not freeze)

25°C	14 days (Unopened)
30°C	5 days (Unopened)
Expiry	15 months after manufacturing date

※ Use LuciPac A3 for Lumitester Smart, PD-30 or PD-20. Do not use it for other models.

## Kikkoman Biochemifa Company

2-1-1 Nishi-Shinbashi, Minato-ku, Tokyo 105-0003 JAPAN  
 TEL : +81-3-5521-5481 FAX : +81-3-5521-5498  
[biochemifa@mail.kikkoman.co.jp](mailto:biochemifa@mail.kikkoman.co.jp)  
<http://www.kikkomana3.com/>  
<http://biochemifa.kikkoman.co.jp/e/>



※ The information contained in this leaflet is subject to change without further notice.  
 © 2018 Kikkoman Corp. (15020181101)

Scan for more info on  
A3 Hygiene Monitoring System

HyServe GmbH & Co KG.  
 Hechenrainerstrasse 24 D- 82449 Uffing am Staffelsee  
 Tel. 0049 – 8846-1344 Fax 0049 – 8846-1342  
[info@hyserve.com](mailto:info@hyserve.com) [www.hyserve.com](http://www.hyserve.com)